

## UMx-CAM Setup Compatibility and Security

### **Q: When do I need to setup the compatibility and security?**

A: When you encounter with following situations, please adjust the setting in **Properties**.


#### ● Not able to activate the Device Control Mode.

It is because UMx-Cam software requires to access to System to control the microscope motor. But system does not allow it.

#### ● When you see “Microscope is Found” but Connect button is white out.

Because the software was build up under Windows XP, therefore, for some new Operation system such as Win 8 and Win 10, you need to change compatibility to Win XP SP3.


#### ● Access Violation:0040E8E5, 07B3571C,

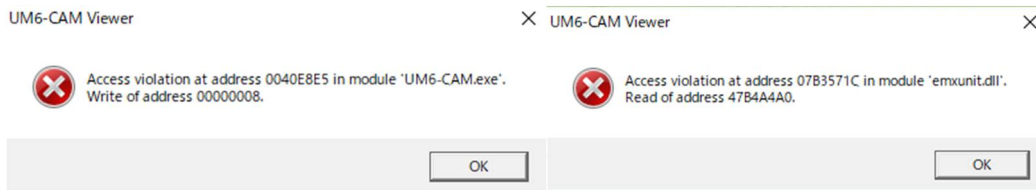


#### ● Unable to turn on UMx-CAM software after Windows OS Update.

Please refer to following steps:

(a) Compatibility : Firstly, Right click the Vitiny software to open **Properties**

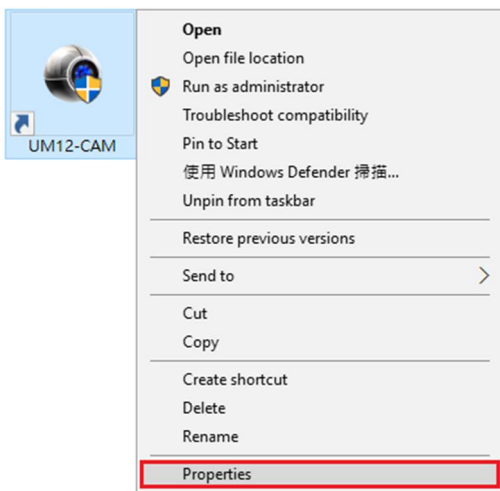

1. To run a program in compatibility mode (See screenshot below)
  - (A) Click on **Compatibility** tab
  - (B) Check the **Run this program in compatibility mode for** box
  - (C) Click on the drop down menu arrow and select Windows XP (Service Pack 3)

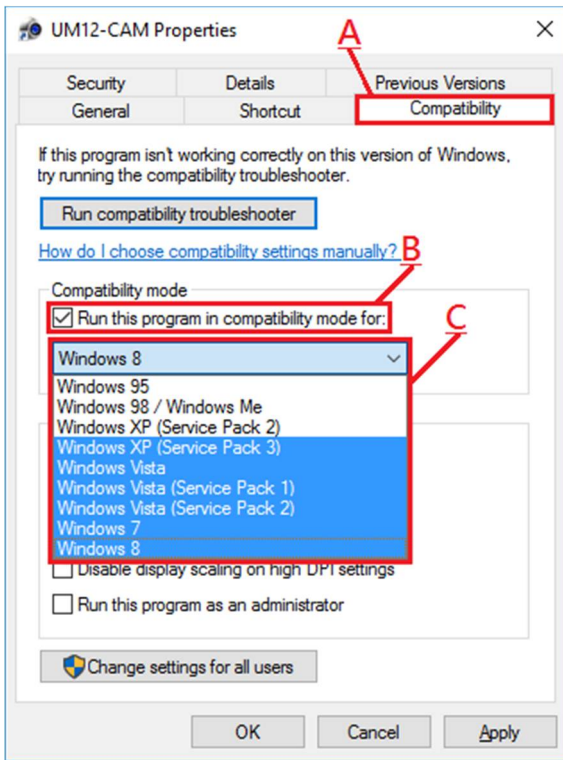


(b) Security: To change the permission for user (See screenshot below)

(D) Click on **Security** tab

(E) Click **Edit**, and (F) choose **\*User (User-PC\Administrator)** to edit permission for user

NOTE: *Name can be User or other according to computer setting.*

(G) Check all the **Allow** column, and click **OK** and **Apply**

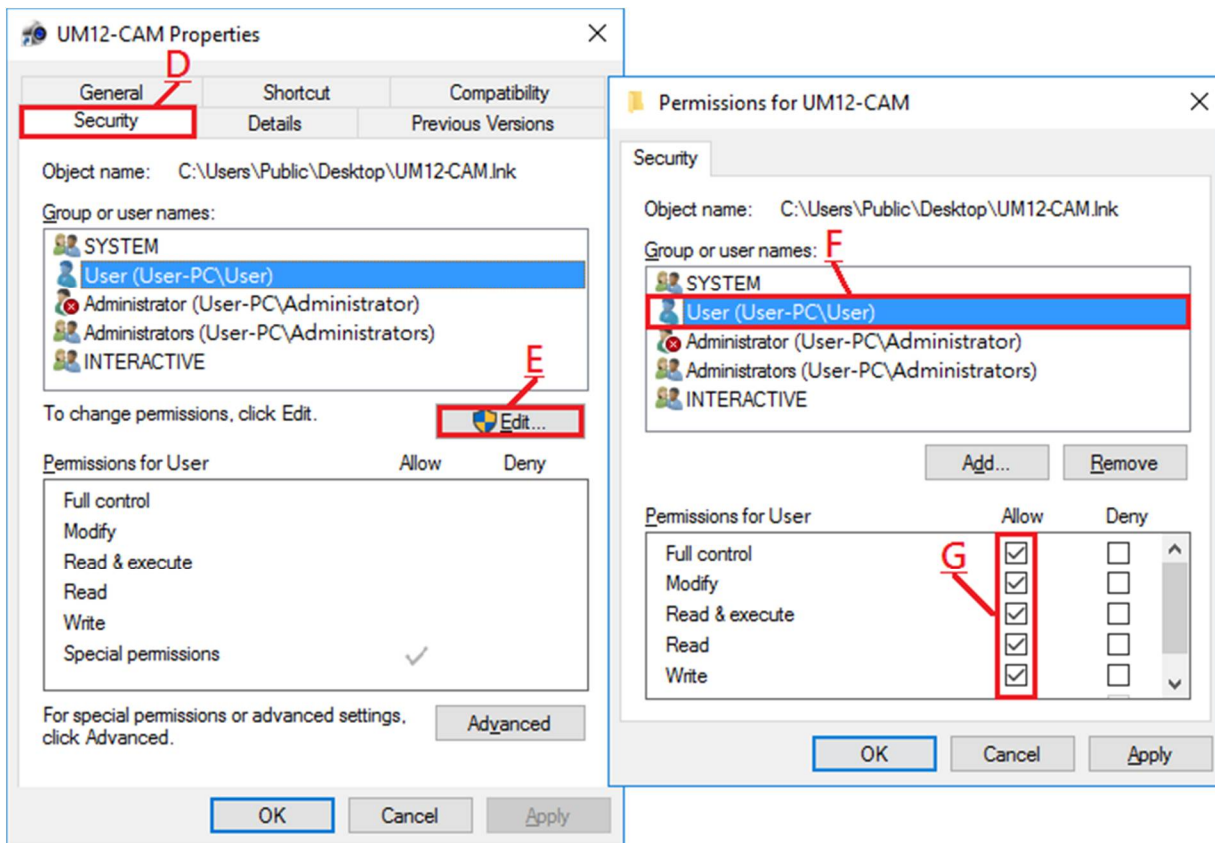

(c) Double check and confirm **OK**.

If you still encounter with connection problem, please contact [mltc@vitiny.com](mailto:mltc@vitiny.com)

