
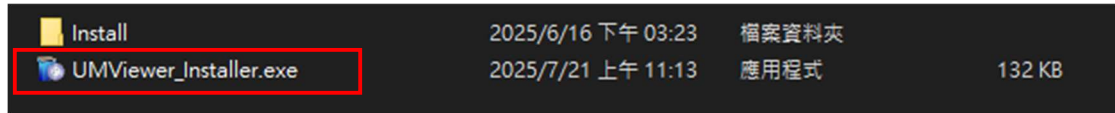
# UMViewer Installation Manual

## Automatic Installation

Select UMViewer\_Installer.exe


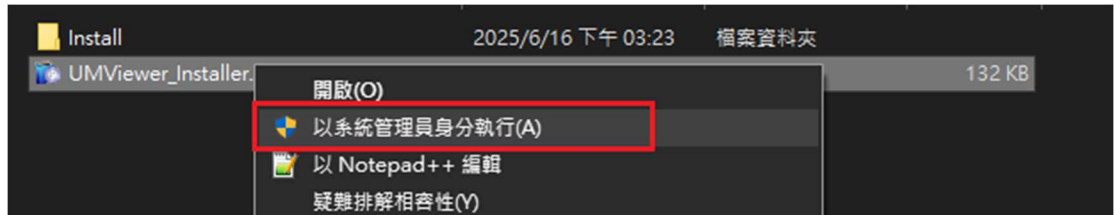
Install UMViewer by Administrator right \_Installer.exe

### A. Click Install Driver

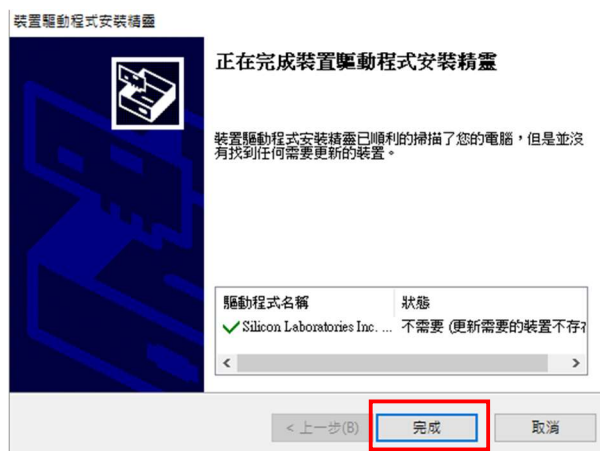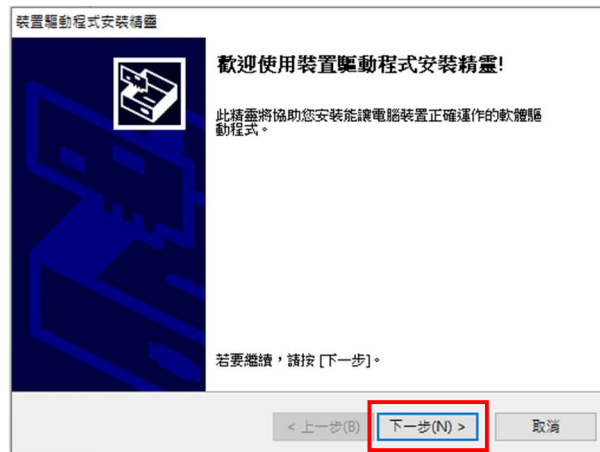If the installation fails, it means the system drive is locked. Please have IT personnel unlock it & try it again.


## B. Click Install Software

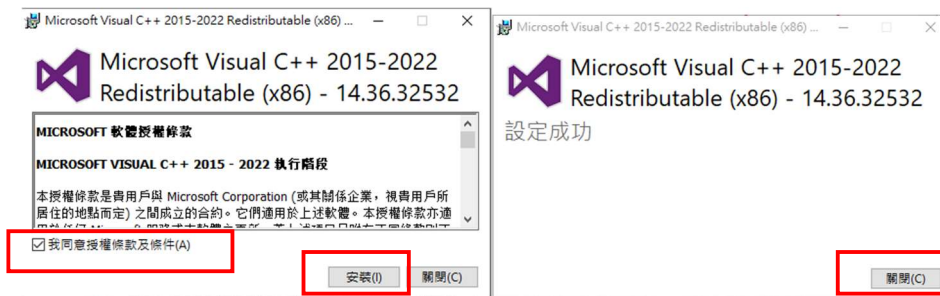
- **Auto detect computer if there is necessary Microsoft Redistributable Package or not. If none, it will remind to install it.**

### 1. Package (1)





Remarks : If the installation fails, please check whether another version is already installed. If so, please close the message and ignore the failure message.

### 2. Package (2)

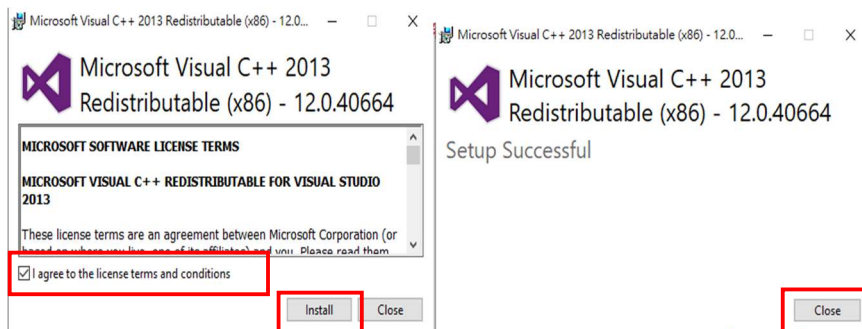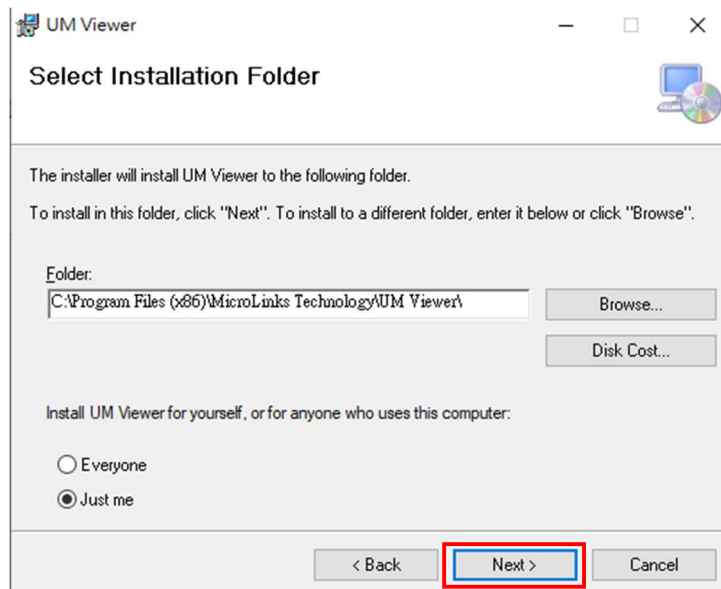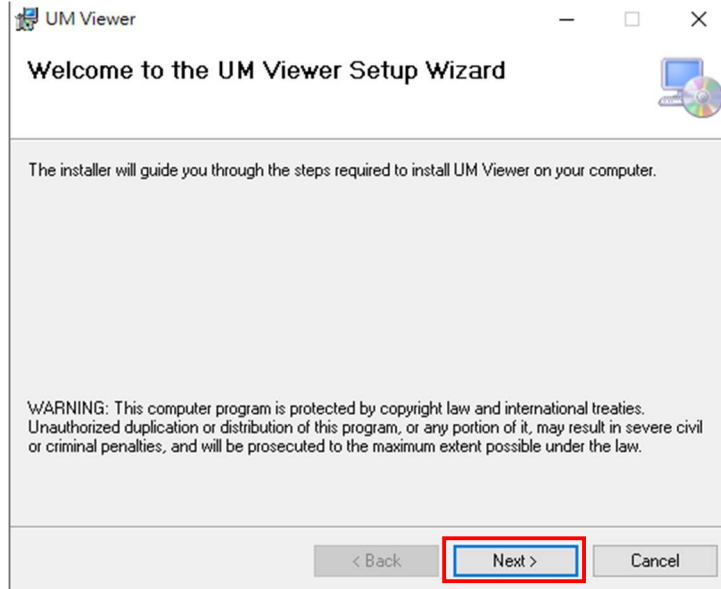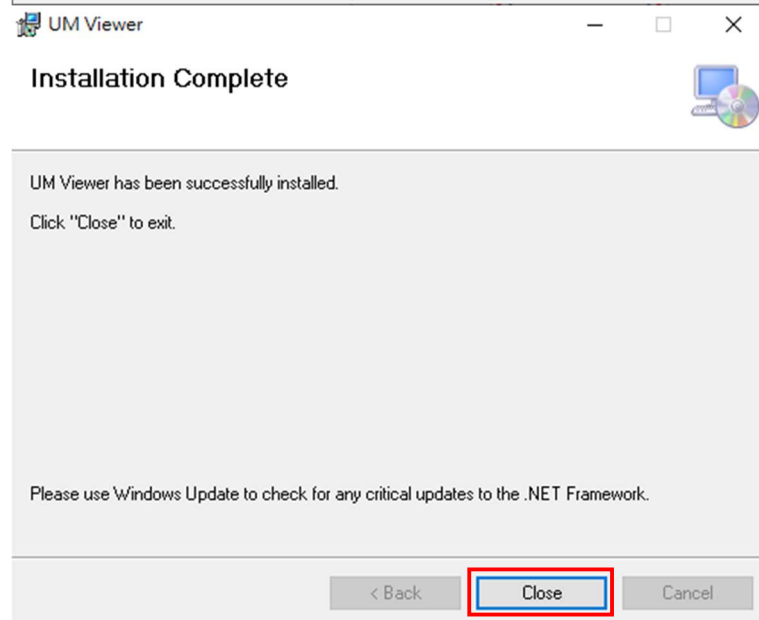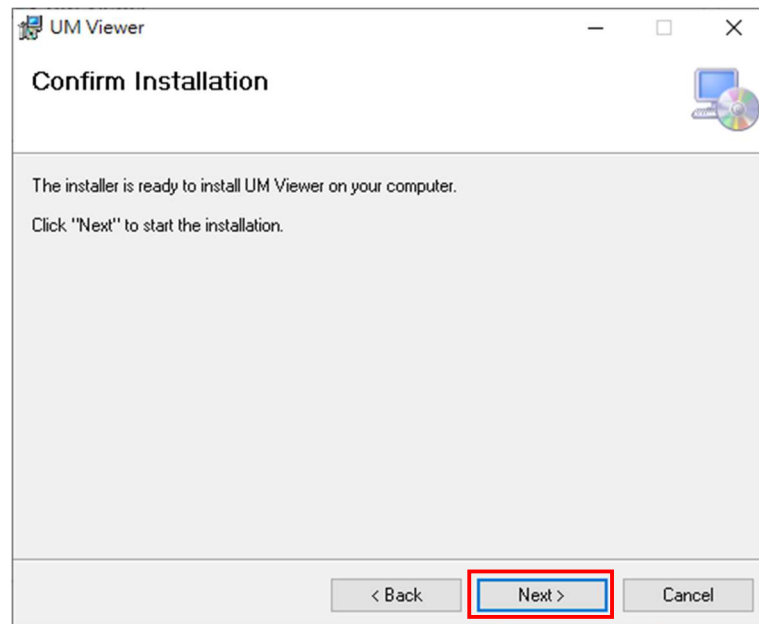Remarks : If the installation fails, please check whether another version is already installed. If so, please close the message and ignore the failure message.

 Installation Finished

## Manual Installation

If automatic installation fails, please use manual installation instead.

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- First, install the driver by clicking the **Driver** folder.

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For **Windows 10** or **Windows 11**, click

10\_

For older operating systems such as **Windows 7** or **Windows 8**, click

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If your operating system is **64-bit**, click

CP210xVCPInstaller\_x64.exe

If your operating system is **32-bit**, click

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

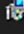
Remarks : If the installation fails, please check whether another version is already installed. If so, please close the message and ignore the failure message.

✚ To install the forwarding **Microsoft Redistributable Package**,  
Click **vcredist\_x86**

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


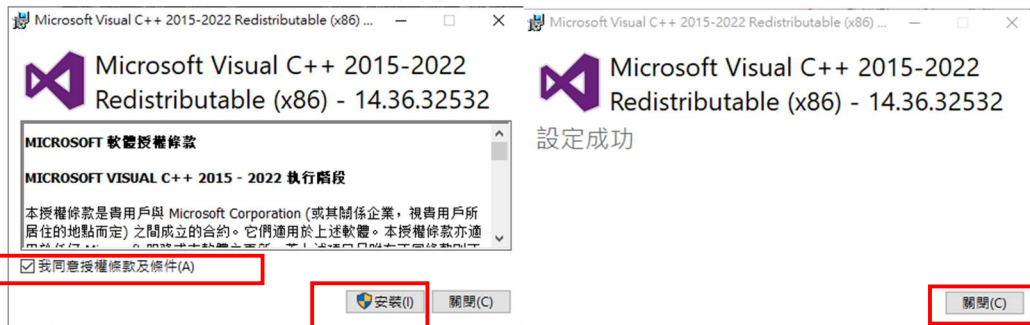
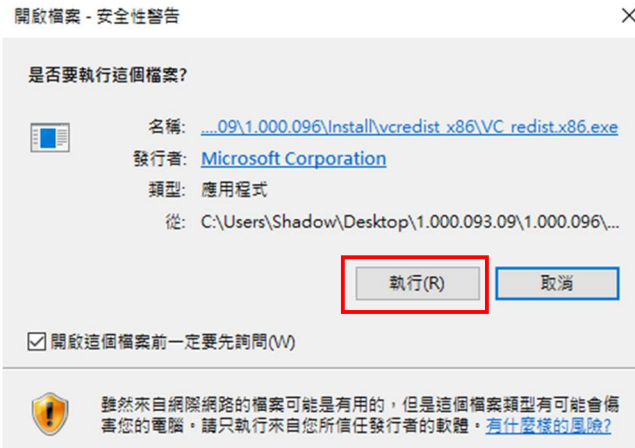
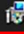
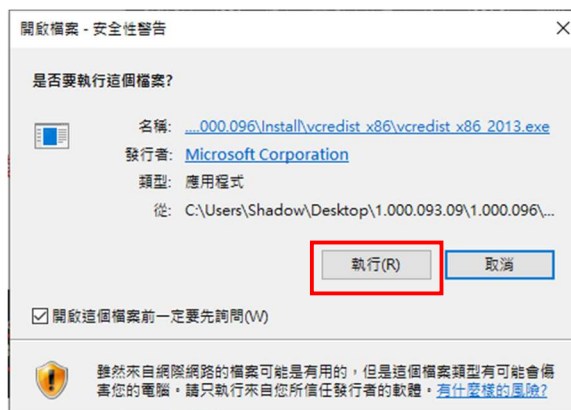
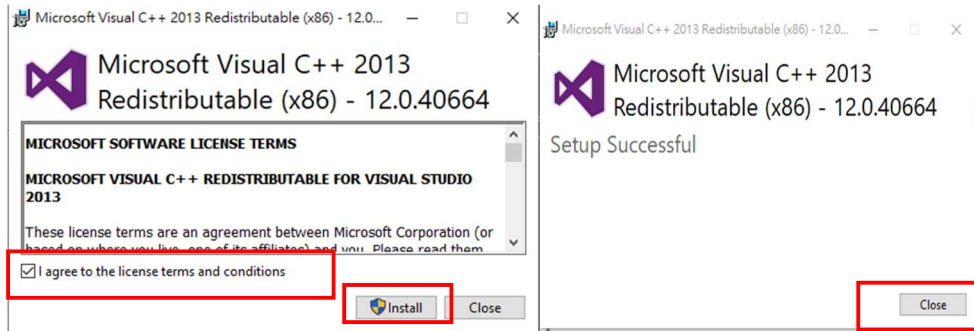
Remarks : If the installation fails, please check whether another version is already installed. If so, please close the message and ignore the failure message.

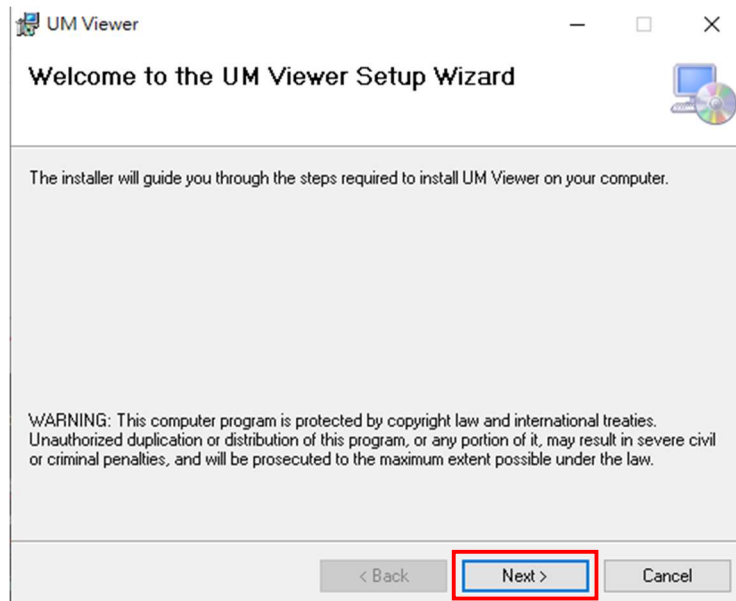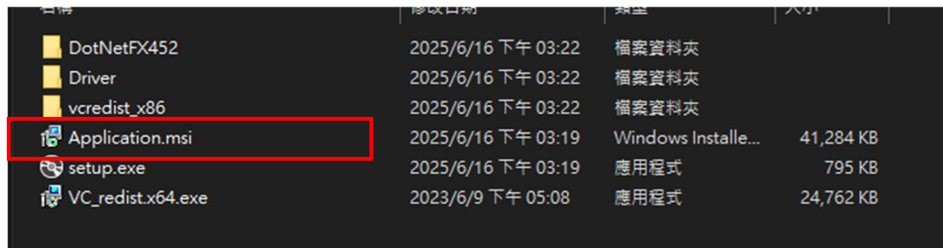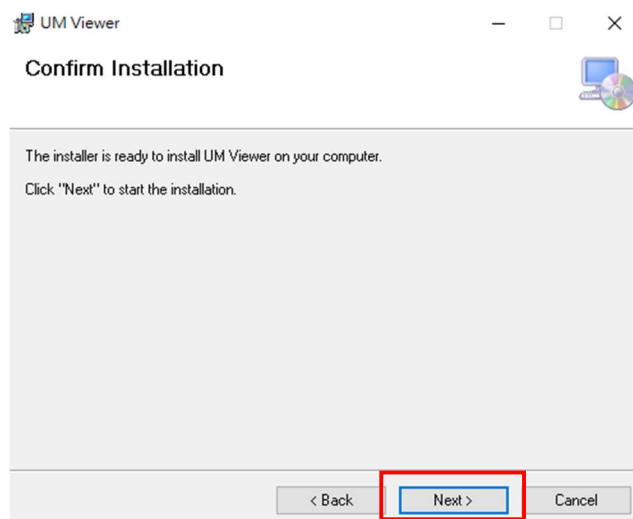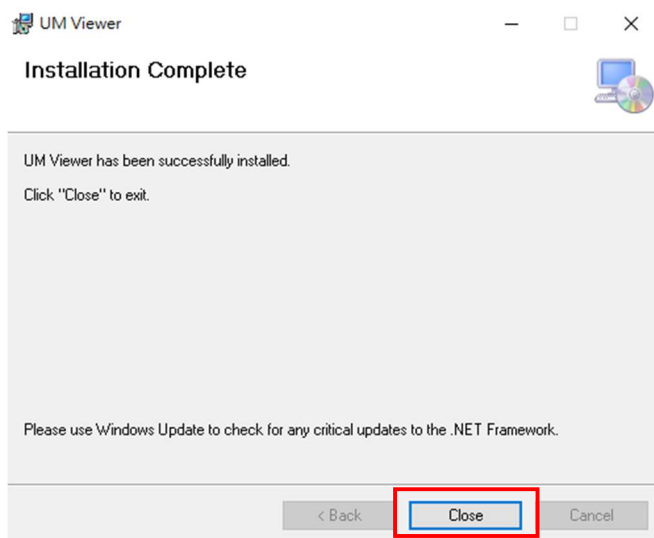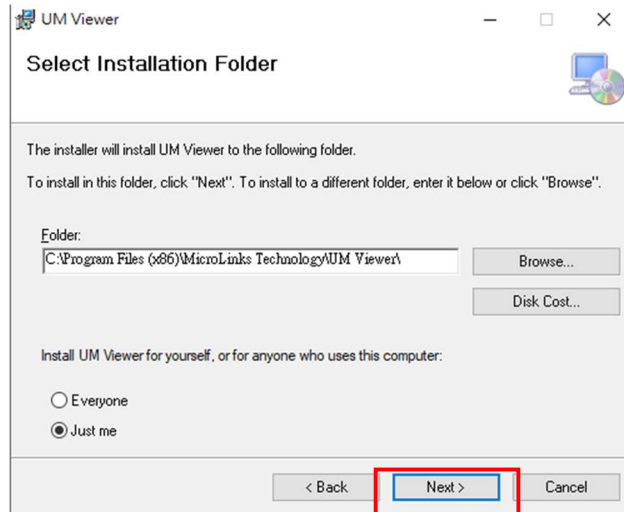To install the UMViewer software, click [Application.msi](#)

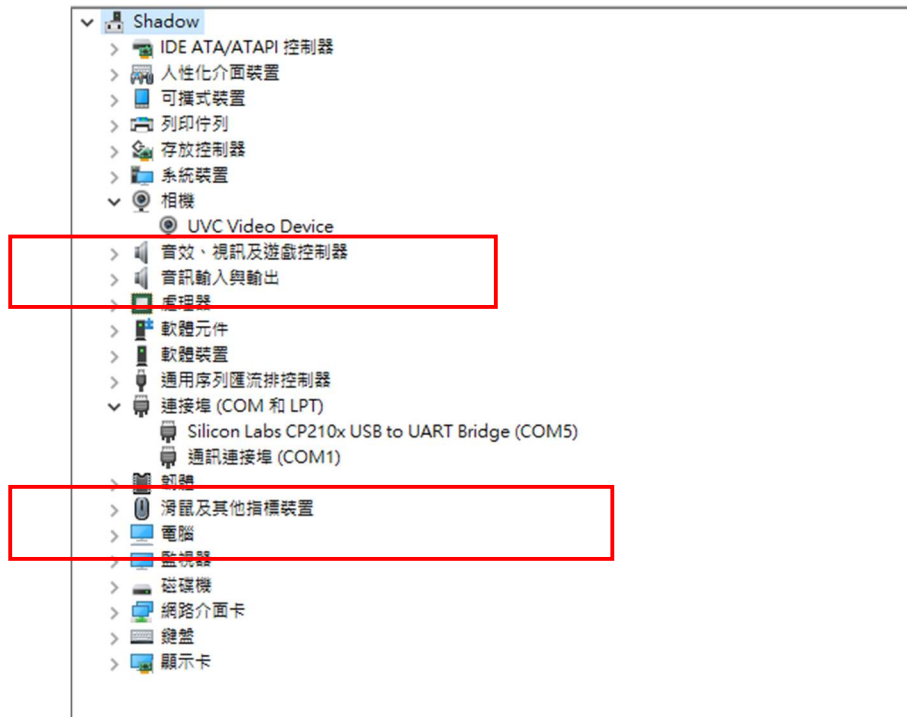
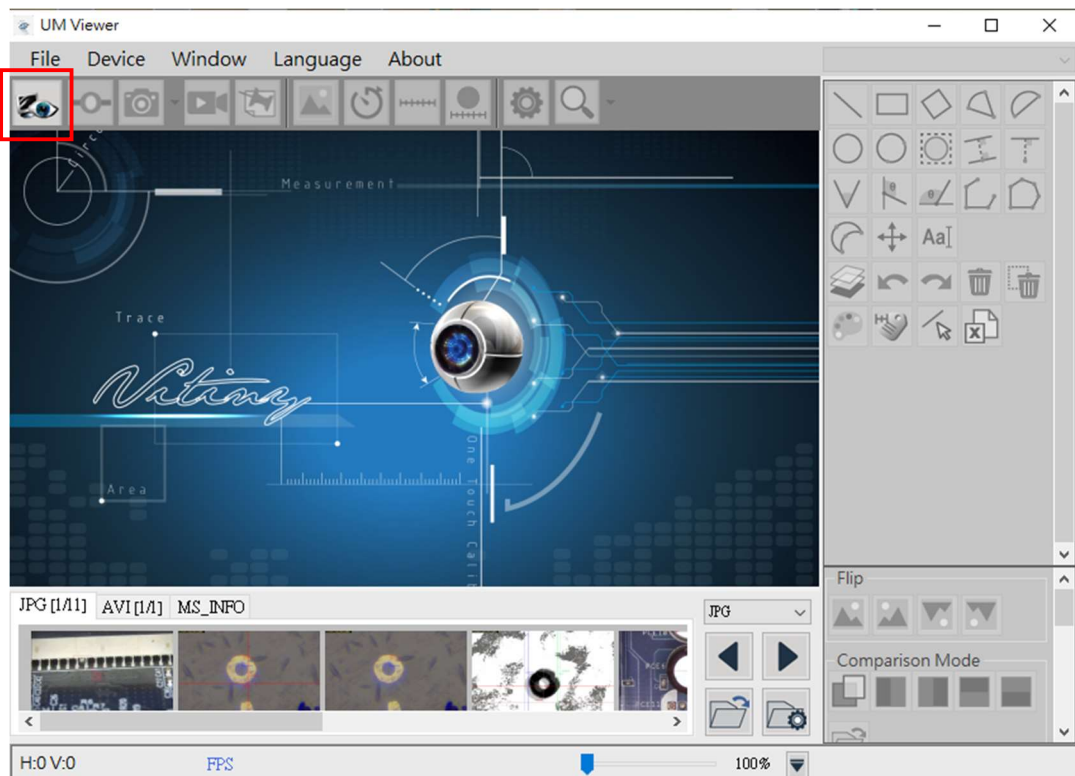
 Installation Completed

# Final check

Connect the microscope > open Device Manager > under Cameras and (COM & LPT), you will see the camera and its driver.

Open the software  this indicates that the device has been detected and can connect.

- ✦ If it is grayed out and cannot connect, click **Device**; the software will perform another search.