
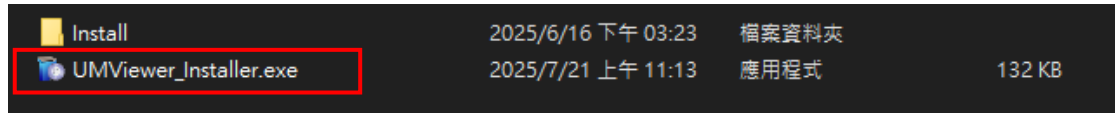
UMViewer Installation Manual

Automatic Installation


Select UMViewer_Installer.exe


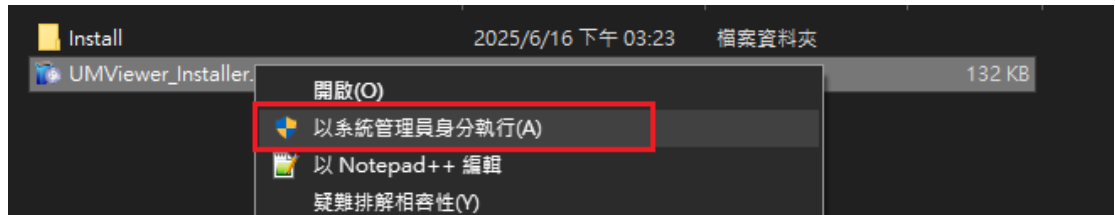
Right click and Install UMViewer by Administrator right _Installer.exe

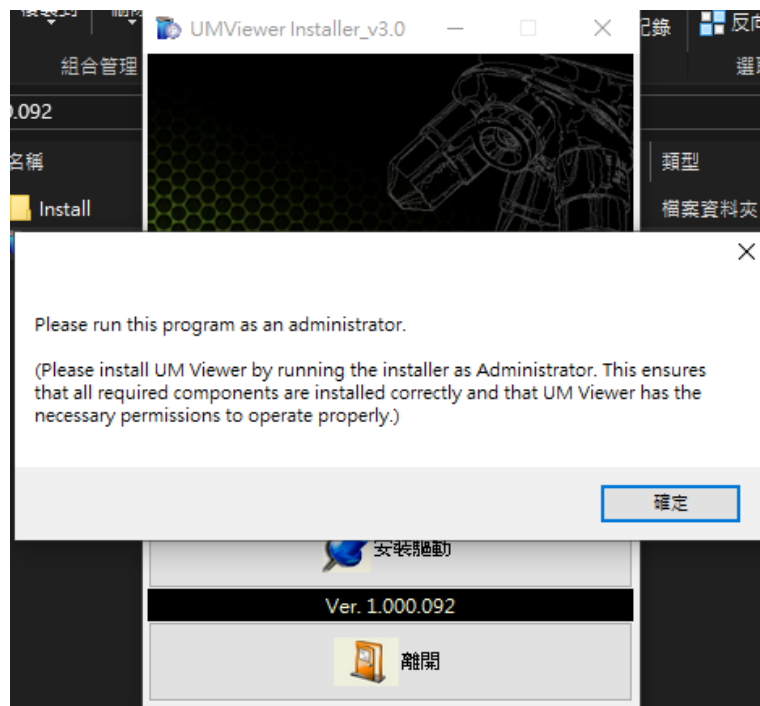
- If it is not selected by administrator right, it will show below message.

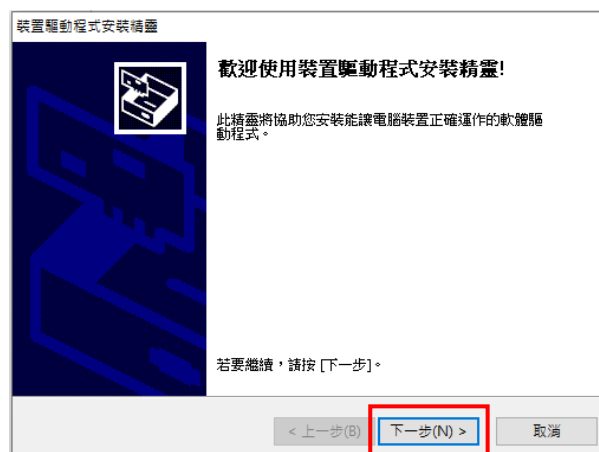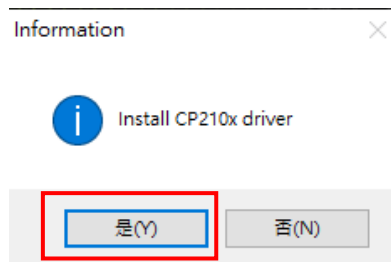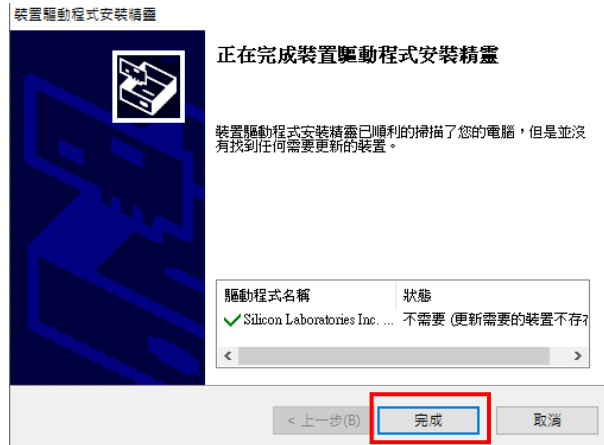A. Click Install Driver


If the installation fails, it means the system drive is locked. Please have IT personnel unlock it & try it again.

B. Click Install Software


- **Auto detect computer if there is necessary Microsoft Redistributable Package or not. If none, it will remind to install it.**

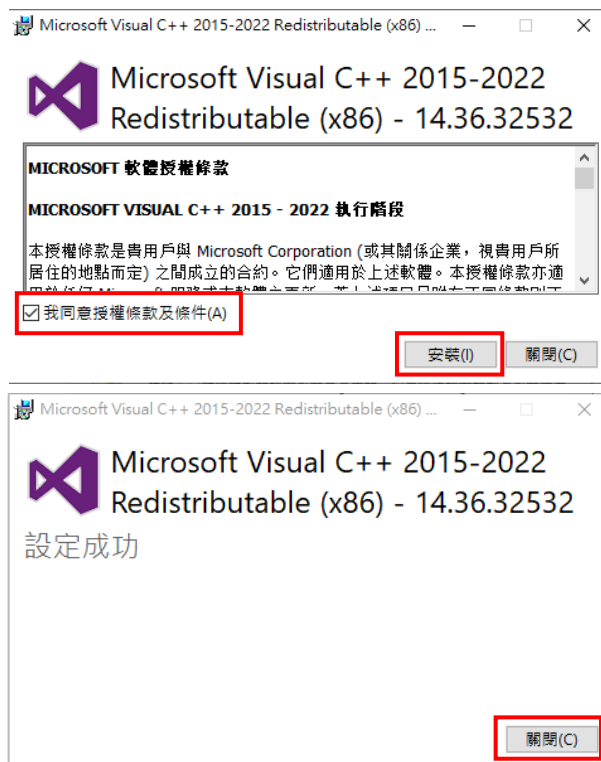
1. Package (1)



Remarks : If the installation fails, please check whether another version is already installed. If so, please close the message and ignore the failure message.

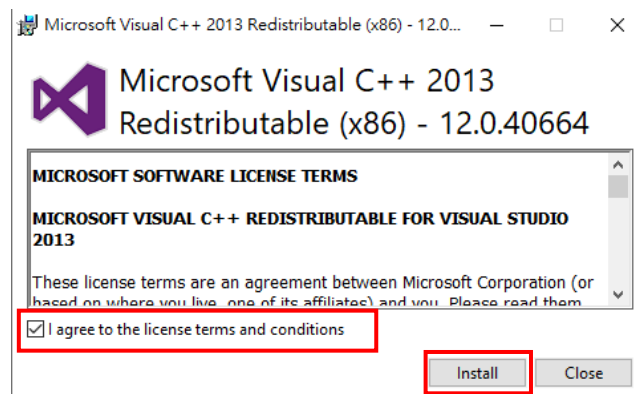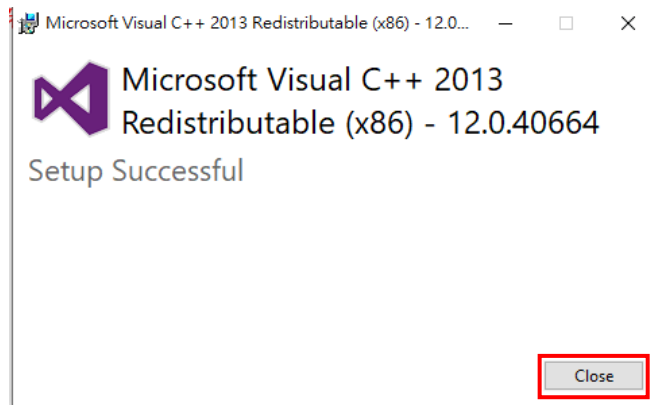
2. Package (2)

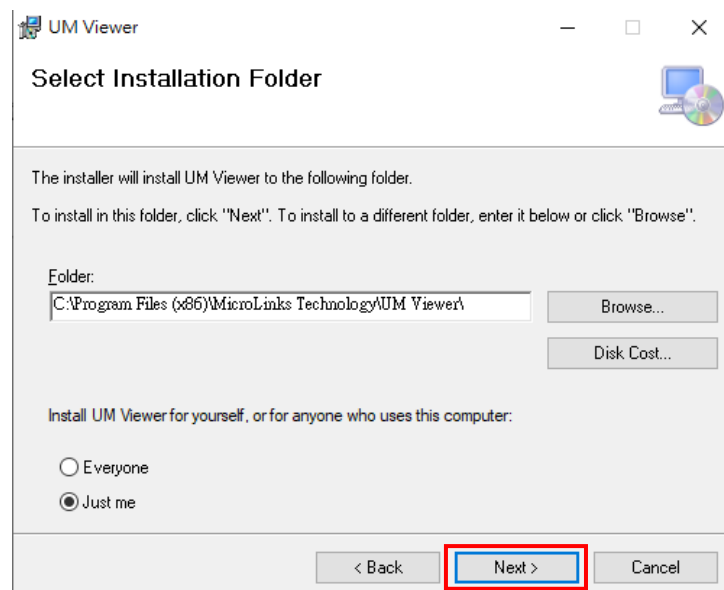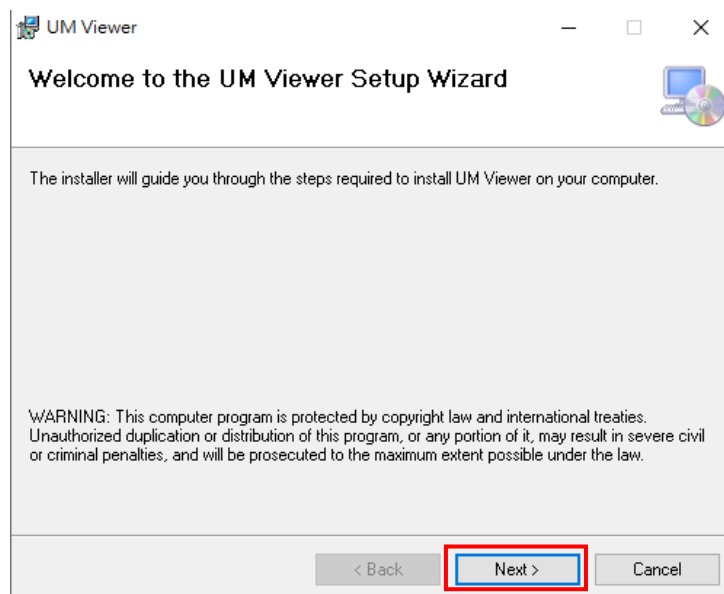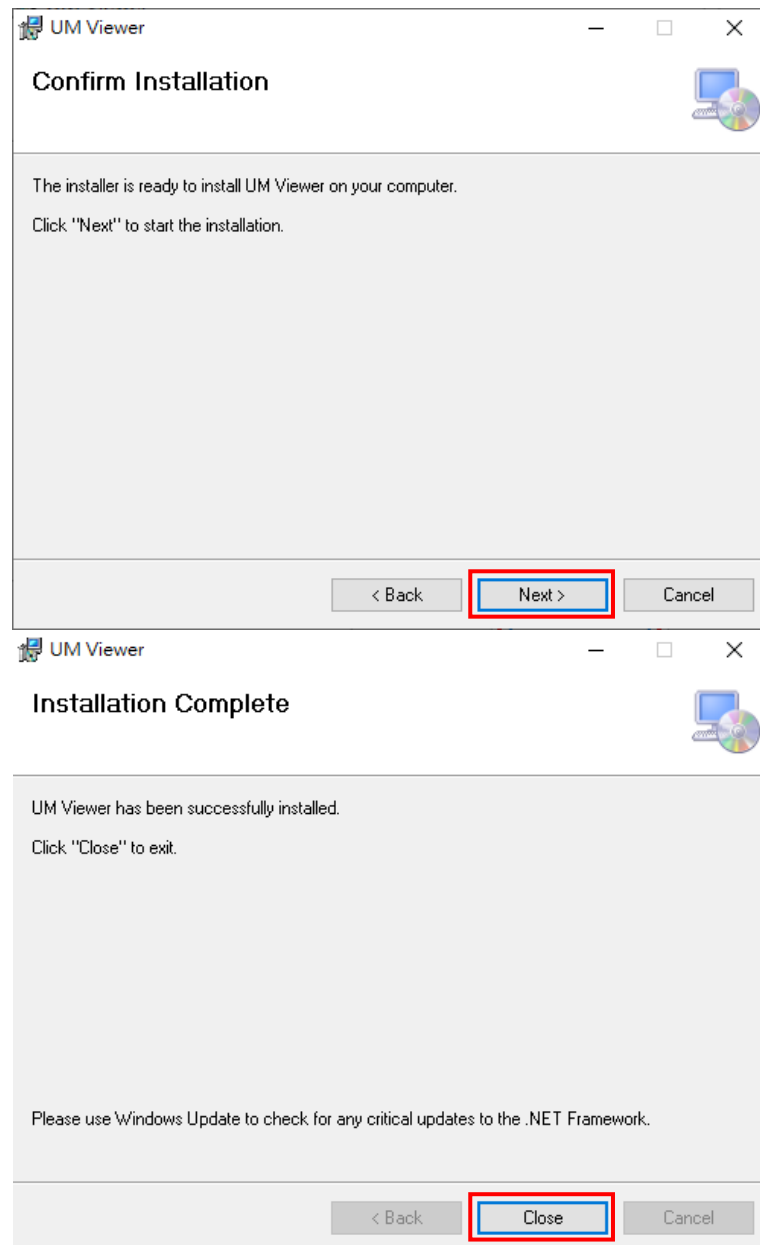Remarks : If the installation fails, please check whether another version is already installed. If so, please close the message and ignore the failure message.

Installation Finished

Manual Installation

If automatic installation fails, please use manual installation instead.

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First, install the driver by clicking the **Driver** folder.

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For **Windows 10** or **Windows 11**, click

 10_

For older operating systems such as **Windows 7** or **Windows 8**, click

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If your operating system is **64-bit**, click

 CP210xVCPInstaller_x64.exe

If your operating system is **32-bit**, click

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
Remarks : If the installation fails, please check whether another version is already installed. If so, please close the message and ignore the failure message.

To install the forwarding package, click **vcredist_x86**

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

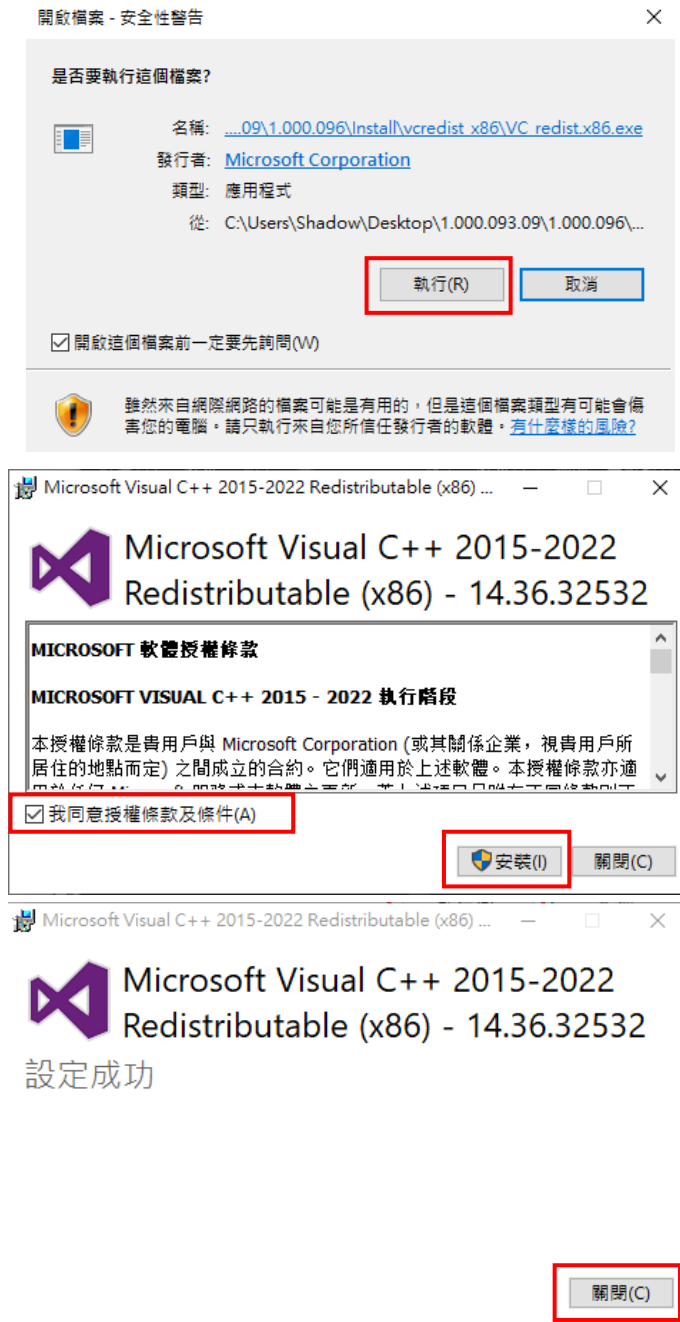
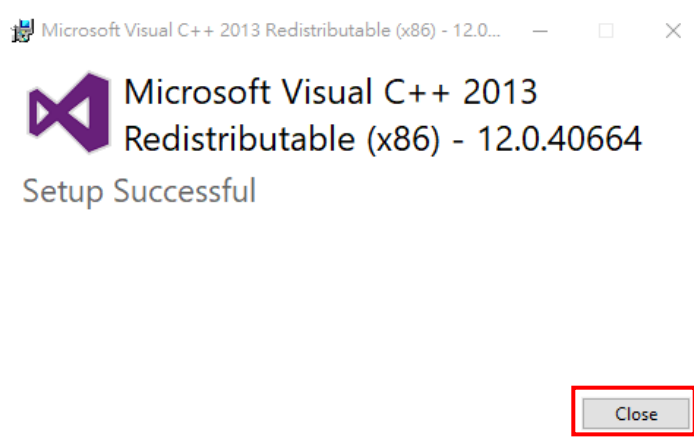
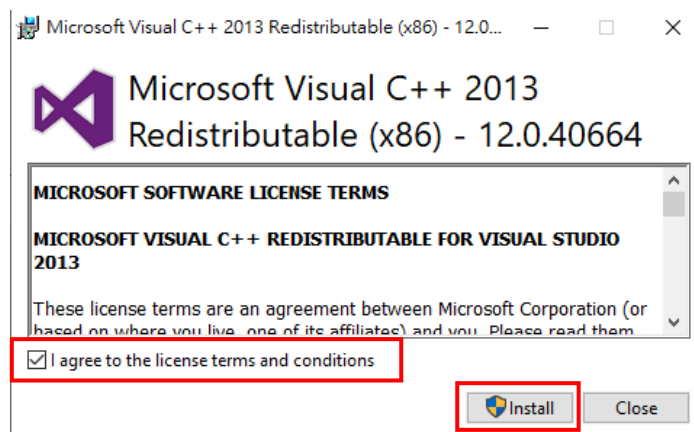
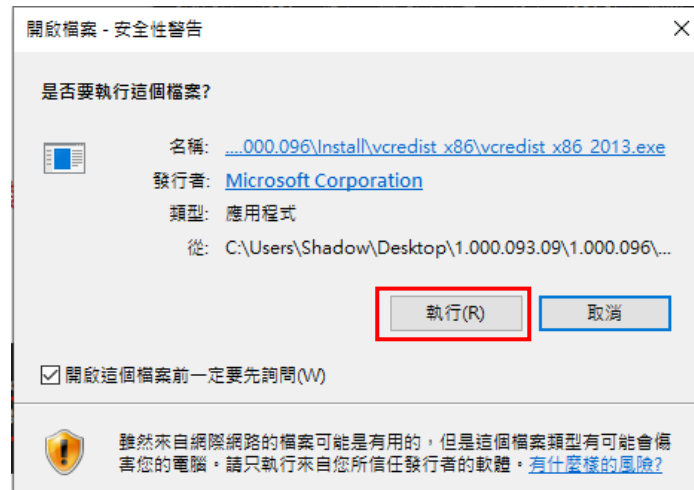
Remarks : If the installation fails, please check whether another version is already installed. If so, please close the message and ignore the failure message.

 To install the UMViewer software, click Application.msi

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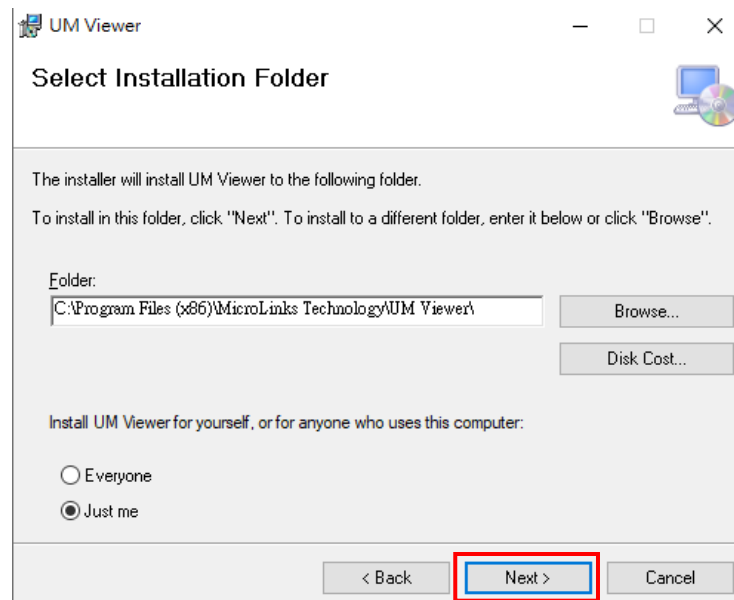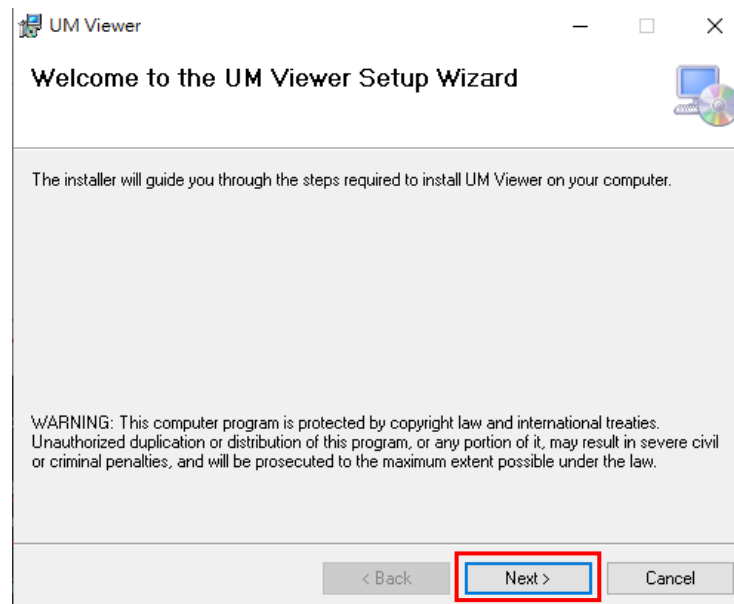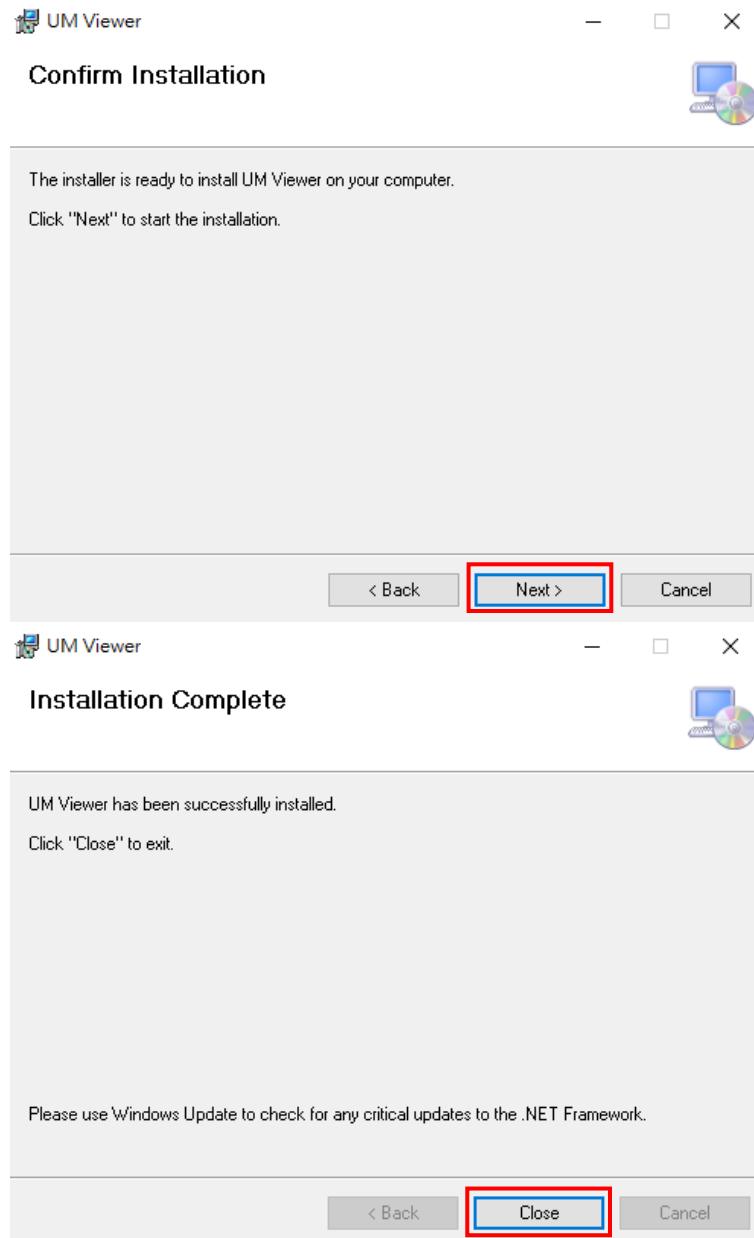
Installation Complete

Final check

Connect the microscope > open Device Manager > under Cameras and (COM & LPT), you will see the camera and its driver.

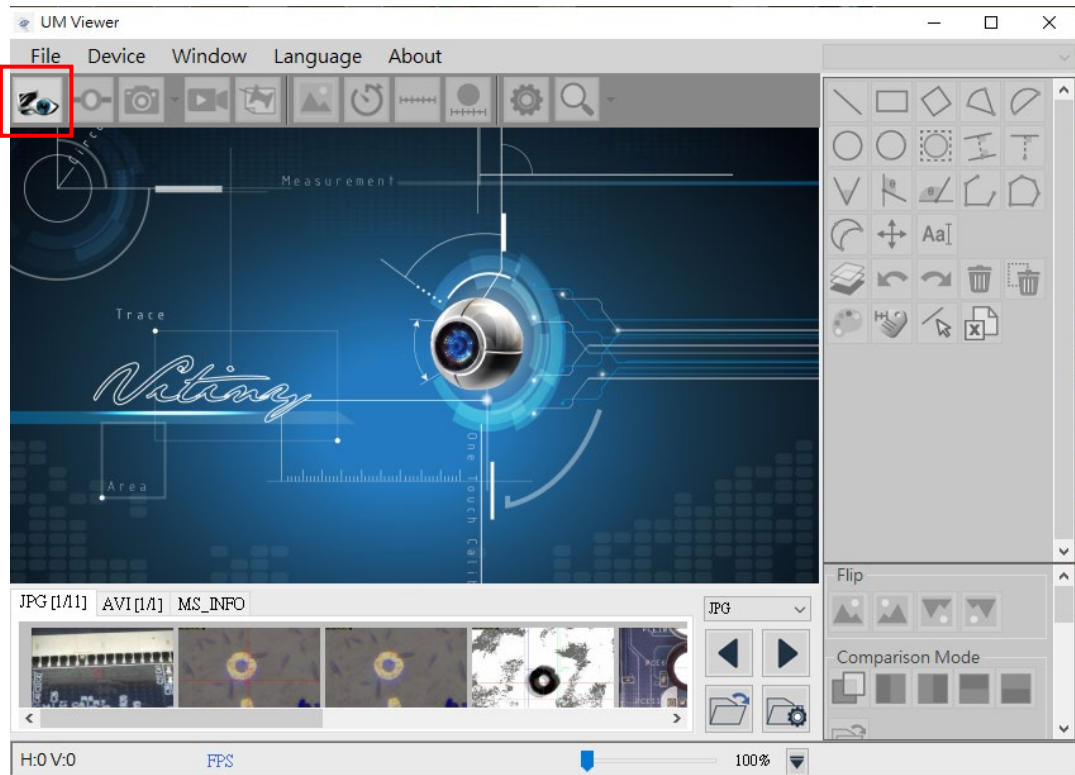Open the software  this indicates that the device has been detected and can connect.

✧ If it is grayed out and cannot connect, click **Device**; the software will perform another search.